



Neutralizes
formaldehyde,
gluteraldehyde
and other aldehyde
solutions!


Neutralin™

neutralizes formaldehyde!

Neutralin™ neutralizes formaldehyde, gluteraldehyde and other aldehyde solutions by converting them into a non-hazardous polymer and water. Fumes and vapors are eliminated and safety is improved. Simply pour used formalin solutions into the five gallon pail containing **Neutralin** and agitate. Neutralization occurs overnight without additional steps or decanting. **Neutralin** is supplied with a Trace Aldehyde Detection Kit, which enables testing of the solution to be sure no residual formaldehyde is left.

Neutralin is prepacked and ready to receive five gallons of 4% formalin solution.

After the **Neutralin** container has been filled and allowed to react overnight, test for the presence of residual formaldehyde using the included **Trace Aldehyde Detection Kit™**. In a small transparent test tube add 0.5ml Aldehyde Test Reagent (from vial provided), 0.5ml Deionized water, and 50µl **Neutralin**

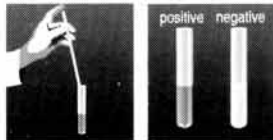

Supernatant (clear liquid from the top of your filled **Neutralin** container). Cap test tube and invert several times. **Negative Test-** Formaldehyde is no longer present and the **Neutralin** reaction complete if the solution remains clear for 5 minutes. **Positive Test-** If formaldehyde is still present, the solution will turn a magenta color **within 5 minutes**. Wait overnight and recheck the solution the next morning. The test should be negative. If the **Neutralin** retests positive (magenta), the solution has been overloaded with formaldehyde. The supernatant (clear liquid) should be decanted into an unused **Neutralin** container. The white solid in the original container should be rinsed with tap water and this water decanted into the new **Neutralin** container. Mix and retest the supernatant for aldehydes. There is sufficient Aldehyde Test Reagent provided to do three tests.

Neutralin System:	Neutralin, 5 gallon pail	\$22.50
Order No. HS-108	including Trace Aldehyde Detection Kit	

305 Patton Drive, Atlanta, GA 30336-1817 (404) 699-2121 FAX: (404) 699-2077

Unit 4, Fleet Business Park, Itlings Lane, Hessle, Hull HU13 9LX, England
0482 646022 FAX: 0482 646013

